

Confocal Start-Up Procedure:

1. Activate anti-vibration system
 - a. (Turn on switch, push “E,” red light on=enabled)
2. Turn on metal halide lamp power supply
3. Turn on PC and microscope switch
4. Log on (password to individual labs = For example, Vsenkalab)
5. Turn on Scanner power switch and wait ~20 seconds before launching software
6. Launch LAS-AF (advanced fluorescence, not LAS-AF hardware)
7. Change objectives using software (has safety- asks about immersion media)
8. If using lasers, turn laser key on at this point and turn on laser switch